
PRODUCT CATALOG 2021 KEOFITT SINGLE-PORT







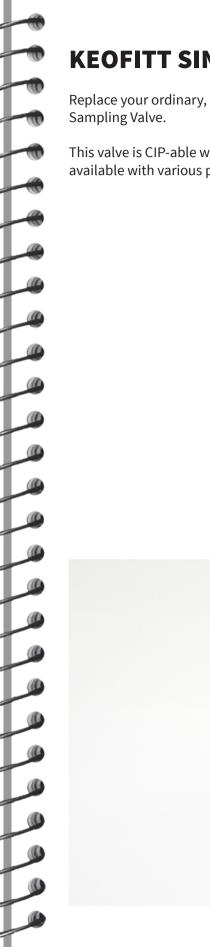

KEOFITT SINGLE-PORT


Replace your ordinary, non-hygienic sampling cocks with the KEOFITT Single-Port

This valve is CIP-able with process CIP and still true to the Keofitt concept, as it is available with various process connections with manual or pneumatic operation.






HOW IT WORKS

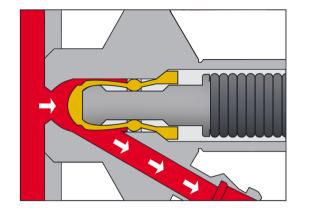

SAMPLING

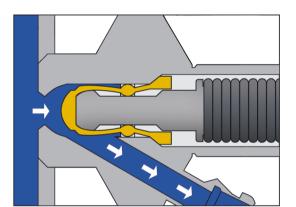

KEOFITT SINGLE-PORT W9[™] SAMPLING VALVE

The KEOFITT Single-Port W9[™] sampling valve is our solution for CIP-able sampling from less demanding applications. If the BASIX[™] does not cover your need due its smaller size, the 9mm outlet on the Single-Port W9 will do the job. With the W9solution you have the added benefit of choosing products from a selected range within the W9 product outline.

This valve is designed for sampling of products with a viscosity of 0-1000 cP with particles up to 3 mm in diameter.

The Single-Port W9 valve is usable for simple, hygienic sampling in all industries for products such as:

Water

Milk


Perfume

Fruit juice

Available in many versions; choose your preferred type of head from our W9 product outline.




KEOFITT SINGLE-PORT W9 - BODIES

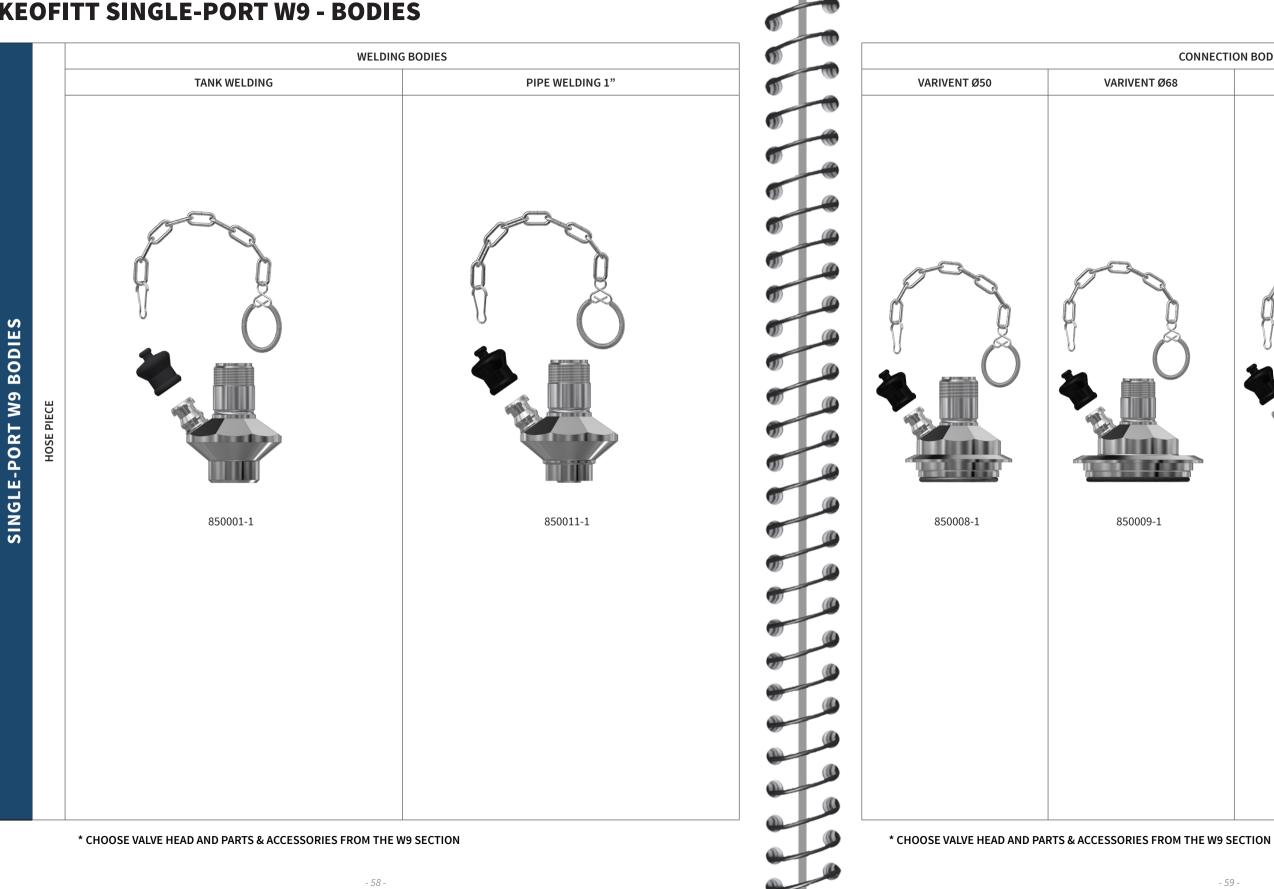


CONNECTION BODIES		
T Ø68	MINI CLAMP	CLAMP 1" & 1½"
9-1	850036-1	850021-1

HOW IT WORKS

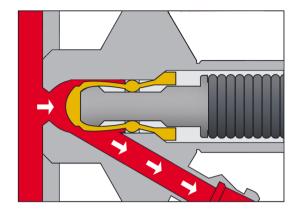

CIP

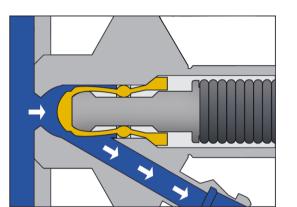

KEOFITT BASIX SAMPLING VALVE

The KEOFITT BASIX[™] Sampling Valve is our basic CIP-able sampling valve for less demanding applications. The KEOFITT BASIX[™] Sampling Valve is our product category for process CIP-able sampling valves for less demanding applications that require no separate inlet for sterilizing or sanitizing media.

This valve is designed for sampling of low viscosity products in the range of 0-100 cP with particles less than 1,5 mm in diameter.

Basix[™] is used for simple, hygienic sampling in all industries for products such as:

Water

Milk

Perfume

Fruit Juice

Available in manual and pneumatic versions this valve is offered complete as a valve body and valve head combination.

Choose from silicone, EPDM or PTFE membranes.



- 60 -

KEOFITT BASIX SAMPLING VALVES MANUAL VALVE PIF TANK WELDING PIPE WELDING MINI CLAMP 1" & 1½" CLAMP TANK WELDING The second BASIX[™] -KEOFITT HOSE PIECE SILICONE 450141 451141 452241 450144 452141 EPDM 450141EPDM 451141EPDM 452141EPDM 452241EPDM 450144EPDM 45 PTFE 450141PTFE 451141PTFE 452141PTFE 452241PTFE 450144PTFE 45

PNEUMATIC VALVE			
IPE WELDING	MINI CLAMP	1" & 1½" CLAMP	
451144	452144	452244	
451144EPDM	452144EPDM	452244EPDM	
451144PTFE	452144PTFE	452244PTFE	

